

PROTOCOL

Protocol for chemotaxis of human peripheral blood B-cell-derived cell lines MAVER-1 (mantle cell lymphoma origin) and MEC-1 (chronic lymphocytic leukemia origin)

CELLS AND REAGENTS: MEC-1 and MAVER-1 cell lines
RPMI1640 medium containing 10 % FBS
Fibronectin
CCL19

Note: All media should be equilibrated before use (overnight at 37 °C in an incubator) to avoid bubble formation during the experiment. This applies to the RPMI1640 medium, Milli-Q water for fibronectin dilution and PBS for cell washes.

Preparation of media

1. Equilibrate RPMI1640 medium containing 10 % FBS overnight at 37 °C in an incubator.
2. The next day, mix part of the medium with the desired chemoattractant (e.g. 200 ng/mL recombinant chemokine CCL19, e.g. 361-MI R&D Systems). To visualize the established gradient, the media can be mixed with a fluorescent marker in 1:20 ratio (e.g. TRITC 20K, 81-001, Gradientech).
3. Fill syringe 1 with control medium (no chemoattractant) and syringe 2 with medium containing the chemoattractant (and fluorescent marker). Keep at 37 °C in the incubator until use.

Pre-coating of the CellDirector 2D surface with fibronectin

Note: Both MEC-1 and MAVER-1 cells are cell lines cultured in cell suspension. MEC-1 cell line tends to adhere weakly to the plate surface, however MAVER-1 do not adhere at all. Therefore, surface coating is crucial for these experiments.

1. Dilute the human plasma-derived fibronectin (e.g. 1918-FN-02M, R&D, stock 1 mg/mL) to a final concentration of 50 µg/mL in Milli-Q water (1:20).
2. Inject 200 µL of the solution to the CellDirector 2D with a pipette. Coat overnight at 37 °C in incubator. Keep the device in a 10 cm petri dish containing a wetted tissue to avoid evaporation of the solution.
3. Next day, replace the fibronectin solution with equilibrated medium - wash twice with 200 µL of the media and then let block the surface for 30 min at 37 °C in the incubator.

Preparation of cells

Note: To simplify the cell-tracking during experiments, the cells can be stained directly before use by membrane-permeable dyes (e.g. Calceins)

1. Harvest the cell suspension from the cell culture flask, spin down (200 x g, 5 min) and wash twice with PBS.
2. *Optional washing step by Calcein Green, AM (C34852, Thermo Fisher Scientific): 10 min of incubation in 2 μ M Calcein Green solution diluted in PBS, followed by 1-2 washing steps. Optimal concentration of the cell suspension: 1-2 x 10⁶ cells/mL.*
3. Resuspend the cells in the equilibrated medium at a final concentration of 1-2 x 10⁶/mL.

Loading the cells

1. Inject slowly the 200 μ L of the cell suspension to the CellDirector.
2. Incubate for 1 hour in the incubator, in a 10 cm Petri dish. Verify proper cell adhesion by microscope.
3. Set the syringes to the pump and place the CellDirector to a holder on your microscope. Carefully attach the tubes with media and mark the tube containing chemoattractant. Attach the tube with vacuum to the device.
4. Set the CO₂ chamber in case of longer experiments.
5. Start the experiment. The injection speed should be optimized. For these cells, initial speed 5 μ L/min was used to set the gradient and then was decreased to 1 μ L/min. In case of low cell adherence, the loose cells can be removed from the device by a transient faster medium flow.
6. Collect images using bright field (or fluorescence) microscopy (10x objective) at 0.5 fpm.

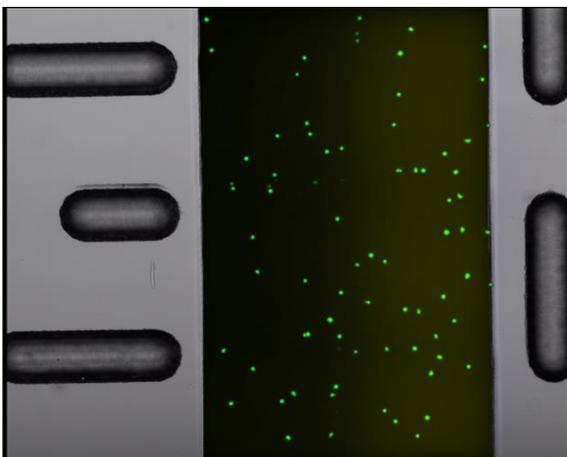


Fig. 1: Calcein-labelled B cell lymphoma-derived cells in CellDirector® 2D gradient channel, with gradient marker.

Protocol modification – switch of treatments (CellDirector® Switch, 71-001, GradienTech)

Note: To change the experimental conditions throughout the running experiment, a CellDirector Switch connected to an additional syringe (e.g. containing inhibitor of the chemotaxis) can be used. Do not forget to assemble this device ahead together with the remaining syringes.

We thank Pavlína Janovská, MSc, and Assoc. Prof. Vítězslav Bryja, PhD, at the Dept. of Experimental Biology, Masarykova Universita, Brno, Czech Republic, for sharing the protocol.